AutoRep™ M Repetitive Pipette

Operating Manual

Please follow step by step!

Mode d'emploi

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RAININ INSTRUMENT, LLC.

E-Mail: pipets@rainin.com

Internet: www.rainin.com

9978 24 · Printed in Germany · 21/1001

RAININ
®

Safety considerations

This instrument may sometimes be used with hazardous materials, operations, and equipment. To address all of the safety problems associated with its use in such applications is beyond the scope of this manual. It is the responsibility of the user of this instrument to consult and establish appropriate safety and health practices and determine the applicability of regulatory limitations prior to use.

1. Observe general instructions for hazard prevention and safety instructions; e.g. wear protective clothing, goggles and gloves. Observe standard laboratory procedures and precautions when pipetting infectious or hazardous samples.
2. Read and follow this Operating Manual.
3. Observe all safety precautions provided by reagent manufacturers. If in doubt, contact the manufacturer or supplier.
4. Use only with samples that do not attack polypropylene and polyethylene.

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5. Never use force on the instrument.
6. Operate only with tip attached (use Encode tips).
7. Avoid touching the tip orifice when working with hazardous samples.
8. Always work in a way which neither endangers the user nor any other person. Dispense samples into a suitable container, avoiding splashing.

Purpose

The AutoRep M is a hand-held dispenser for easy repetitive dispensing of samples. After selecting an Encode tip from a range of sizes and setting the stroke length, volumes from 10 μL to 5 mL can be dispensed with the highest precision and accuracy within the following limitations.

- 15 °C to 40 °C (59 °F to 104 °F)
(instrument and reagents
- other temperatures on request)
- vapor pressure up to 500 mbar
- viscosity:
 - 20 mPa s with 50 mL Encode tips
 - 260 mPa s with 5 mL Encode tips
 - 977 mPa s with 1.25 mL Encode tips

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Operating exclusions

During proper operation, the dispensed sample only comes in contact with the tip and not with the AutoRep M.

Never dispense samples which attack polypropylene or polyethylene.

Do not use aggressive cleaning agents to clean the pipette.

Note:

The instrument and the Encode tips are not autoclavable (121 °C / 250 °F). See sterile Encode tips.

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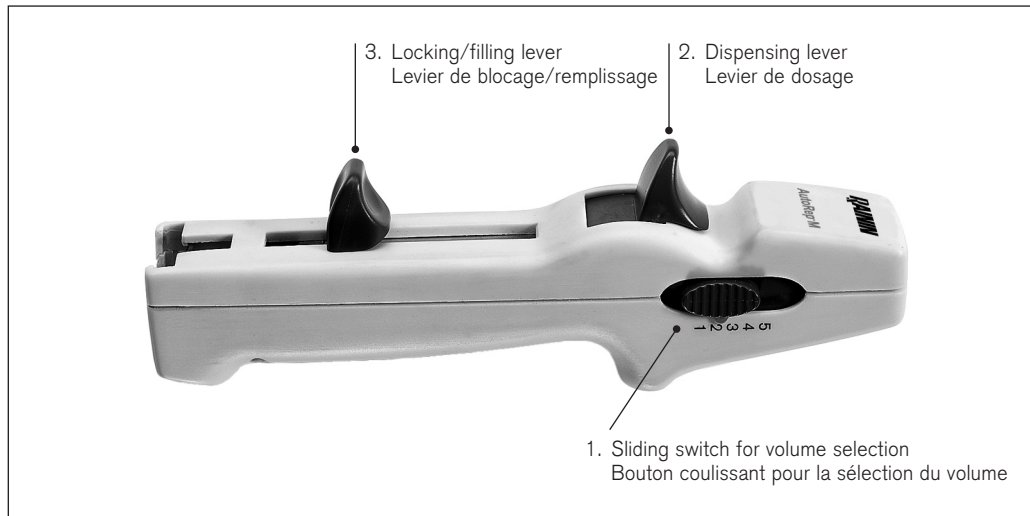
Operating elements

1. Sliding switch for volume selection
2. Dispensing lever
3. Locking/filling lever

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
Operation

Mounting the tip

1. Push the locking/filling lever to the lower stop.
2. Pull out the locking/filling lever slightly and swing it out.
3. Insert the Encode tip.
4. Swing the locking/filling lever back to lock the tip.

Note:

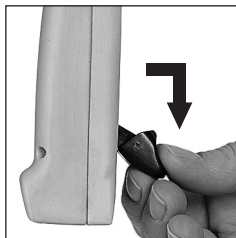To mount 25 mL and 50 mL Encode tips, use the adapter supplied with the tips. The reusable adapter twists on and off.

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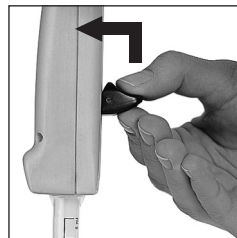
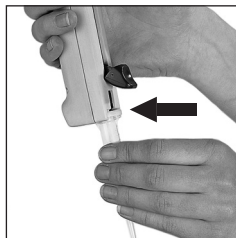
Adjusting the desired volume

Dispense volumes for the AutoRep M are shown in Table 1 below. The table is a matrix of all sizes of Encode tips and all stroke settings.

The dispensed volume is obtained by multiplying the smallest volume unit (printed on the Encode tip) by the stroke setting.

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Filling the Encode tip

1. Immerse the orifice of the Encode tip to a depth of 3 - 10 mm into the sample.
2. Raise the locking/filling lever slowly to prevent cavitation.

Discard the first dispensing step.

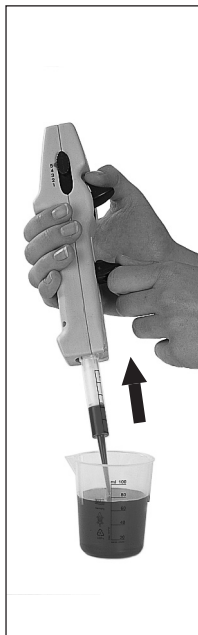
Any small air bubble near the piston can be ignored since a locking mechanism prevents the residual stroke from being dispensed.

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Repetitive dispensing

1. Verify the volume setting.
Wipe off excess sample on the Encode tip around the orifice with a lint-free tissue.
2. Place the orifice of the Encode tip against the inner wall of the vessel.
3. Dispense the sample by pressing down the dispensing lever completely each time.
4. Take care to dispense smoothly and evenly.

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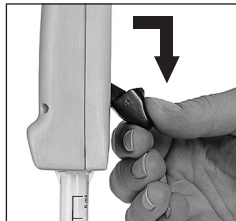
Removing the Encode tip

1. Hold the AutoRep M over a vessel.
2. Empty the Encode tip by pushing the locking/filling lever to its lower stop.
3. Pull out the locking/filling lever slightly and swing it out.
4. Remove the Encode tip.

The adapter required for 25 mL and 50 mL Encode tip is reusable.

Note:

The tightness of the seal of the Encode tips is not guaranteed for repeated use with high viscosity media.

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Troubleshooting

Trouble	Possible cause	Corrective action
Encode tip cannot be inserted	Locking/filling lever is not pushed all the way down and is not swung out The Encode tip piston is not fully inserted	Push the locking/filling lever to its lower stop and swing it out Push the piston of the Encode tip fully into the cylinder.
Locking/filling lever cannot be pushed up	Locking/filling lever is not fully pushed in	Make sure locking/filling lever is fully pushed in
AutoRep M dispenses unevenly	Sliding switch is not locked into position	Lock sliding switch into position
Encode tip drips	Encode tip seal worn Tip loosely mounted	Replace Encode tip Push tip on firmly

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Maintenance

The AutoRep M is calibrated at the factory, and is maintenance-free. It may be cleaned with water or isopropyl alcohol.

The AutoRep M contains no user-servicable parts. Never attempt to disassemble the AutoRep M.

Technical data

Measuring results were obtained by using original Encode tips from Rainin. According to ISO 8655, values are based on the maximum volumes for each tip size. Sample is distilled water at 20 °C / 68 °F.

  20°C
EX

Volume	Accuracy	Coefficient of variation
Encode tip 0.5 mL (10 - 50 µL)	± 0.8%	≤ 1.4% to ≤ 0.6%
Encode tip 1.25 mL (25 - 125 µL)	± 0.8%	≤ 0.8% to ≤ 0.3%
Encode tip 2.5 mL (50 - 250 µL)	± 0.7%	≤ 0.8% to ≤ 0.2%
Encode tip 5 mL (100 - 500 µL)	± 0.5%	≤ 0.6% to ≤ 0.2%
Encode tip 12.5 mL (250 - 1250 µL)	± 0.3%	≤ 0.4% to ≤ 0.2%
Encode tip 25 mL (500 - 2500 µL)	± 0.3%	≤ 0.4% to ≤ 0.2%
Encode tip 50 mL (1000 - 5000 µL)	± 0.3%	≤ 0.4% to ≤ 0.15%

The AutoRep M can also be used with other compatible tips.

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  20°C
EX

Volume	Exactitude	Coefficient de variation
Pointe Encode tip 0.5 mL (10 - 50 µL)	± 0.8%	≤ 1.4% to ≤ 0.6%
Pointe Encode tip 1.25 mL (25 - 125 µL)	± 0.8%	≤ 0.8% to ≤ 0.3%
Pointe Encode tip 2.5 mL (50 - 250 µL)	± 0.7%	≤ 0.8% to ≤ 0.2%
Pointe Encode tip 5 mL (100 - 500 µL)	± 0.5%	≤ 0.6% to ≤ 0.2%
Pointe Encode tip 12.5 mL (250 - 1250 µL)	± 0.3%	≤ 0.4% to ≤ 0.2%
Pointe Encode tip 25 mL (500 - 2500 µL)	± 0.3%	≤ 0.4% to ≤ 0.2%
Pointe Encode tip 50 mL (1000 - 5000 µL)	± 0.3%	≤ 0.4% to ≤ 0.15%

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Return for repair

Important:

For safety reasons, only clean and decontaminated instruments will be accepted for inspection or repair.

- Thoroughly clean and decontaminate the instrument.
- Call 800-543-4030 for a Return Authorization Number
- Send the AutoRep M in a suitable packing container to Rainin Service, Rainin Road, Woburn, MA 01801, describing:
 - The nature of the problem
 - What samples have been used

Note:

No warranty claims can be considered if non-authorized repairs (repairs other than by trained Rainin technicians) have been made or attempted.

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To Order, Call 800-472-4646

Manual repetitive pipette
with wall support

Wall support

Cat. No.: AR-M2

Cat. No.: AR-HU

Encode Positive-displacement tips

Capacity (mL)	package of	Cat. No.: (non sterile)	Cat. No.: (sterile**)
0.5 mL	100	ENC-500	ENC-500S
1.25 mL	100	ENC-1250	ENC-1250S
2.5 mL	100	ENC-2500	ENC-2500S
5 mL	100	ENC-5ML	ENC-5MLS
12.5 mL	100	ENC-12ML	ENC-12 MLS
25 mL*	50/25***	ENC-25ML	ENC-25MLS
50 mL*	25	ENC-50ML	ENC-50MLS

Encode Positive-displacement tips

assorted, 20 each:
0.5, 1.25, 2.5, 5.0 and 12.5 mL

Cat. No.: ENC-SET

Adapter for 25 and 50 mL Encode tips, PP, autoclavable

package of	Cat. No.: (non-sterile)	package of	Cat. No.: (sterile**)
10	ENC-10ADP	1	ENC-ADPS

* incl. 1 adapter, ** sterile/free of endotoxin – single wrapped,
*** Encode tips 25 mL: non sterile pack of 50 / sterile pack of 25

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